## **Freezer Temperature Log**

Month/Year	(Days 1-15	)

Freezer Location \_\_\_\_\_

## Record CURRENT, MIN, and MAX temperatures twice a day. Complete steps 1-3

Step 1

C0

Write your initials and time of day.

Provider Name:	 
Pin number:	

File this log at the end of the month and keep it for 3 years

Staff Initials	Example																															
Day of Month	1		1	4	2		3	2	ł		5	l	5	7		8		9		10		11		12		13		14		15		
	7:30																															
Time	am	am	pm																													
Current	-3																															
MIN	-9																															
MAX	1																															



Step 3

Read the current temperature on thermometer display and record on the temp log. Review the minimum and maximum temperatures and record on the temp log.

Evaluate temperatures. If temperatures are in the DANGER zone (see zone below),



Review/Mark Button: Press to see min/max temperatures

Start/Clear/Stop Button: Press to return to normal mode

-10° & lower -9° -8° -7° -6° -5° -4° -3° -2° -1° 0° 1° 2° 3° 4° 5°	6° 7° 8° 9° 10° 11° 12° 13° & higher
THESE TEMPERATURES ARE OK	DANGER ZONE - TOO HOT
<ul> <li>Vaccine Storage Best Practices</li> <li>Store vaccine at 0° or lower to reduce chance of temperatures exceeding the recommended range</li> <li>Place thermometer in a central area of the unit directly with the vaccines</li> <li>Do not place thermometer or vaccine in the door, near or against the walls, close to vents, or on the floor of the unit</li> </ul>	<ul> <li>If you ever see temps in this danger zone:         <ul> <li>Alert your supervisor immediately</li> <li>Wait no more than 60 min and check to see if the temperature has returned to the proper range</li> <li>If temps are still in DANGER ZONE after 60 minutes, download the data and email to chicagovfc@cityofchicago.org</li> <li>Document the date and actions you take on the Vaccine Storage Troubleshooting Record</li> </ul> </li> </ul>

take immediate action described below.



## **Freezer Temperature Log**

Month/Year	(Days 16-31)	1
------------	--------------	---

Refrigerator Location \_\_\_\_\_

## Provider Name:

Step 1

Write your initials and time of day.

Pin number			
	File this log at the end of the month and keep	it for 3	years

Staff	Example RC																																
Day of Month	15	1	.6	1	7	1	8	1	9	2	0	2	1	2	22	2	3	2	24	2	25	2	6	2	7	2	8	2	9	3	0	3	1
Time	7:30																																
	am	am	pm																														
Current	-3																																
MIN	-9																																
MAX	1																																



Read the current temperature on thermometer display and record on the temp log. Review the minimum and maximum temperatures and record on the temp log.

Record CURRENT, MIN, and MAX temperatures twice a day. Complete steps 1-3

Step 3

Evaluate temperatures. If temperatures are in the DANGER zone (see zone below), take immediate action described below.



Review/Mark Button: Press to see min/max temperatures

Start/Clear/Stop Button: Press to return to normal mode

-10° & lower -9° -8° -7° -6° -5° -4° -3° -2° -1° 0° 1° 2° 3° 4° 5°	6°7°8°9°10°11°12°13°& higher
THESE TEMPERATURES ARE OK	DANGER ZONE - TOO HOT
<ul> <li>Vaccine Storage Best Practices</li> <li>Store vaccine at 0° or lower to reduce chance of temperatures exceeding the recommended range</li> <li>Place thermometer in a central area of the unit directly with the vaccines</li> <li>Do not place thermometer or vaccine in the door, near or against the walls, close to vents, or on the floor of the unit</li> </ul>	<ul> <li>If you ever see temps in this danger zone:         <ul> <li>Alert your supervisor immediately</li> <li>Wait no more than 60 min and check to see if the temperature has returned to the proper range</li> <li>If temps are still in DANGER ZONE after 60 minutes, download the data and email to chicagovfc@cityofchicago.org</li> <li>Document the date and actions you take on the Vaccine Storage Troubleshooting Record</li> </ul> </li> </ul>